# Characterize your reaction at a glance. RTCal a new real time calorimetry technique

Fabio Visentin, Olivier Ubrich, Mettler-Toledo GmbH, AutoChem, Sonnenbergstrasse 74, P.O. Box, CH-8603 Schwerzenbach, Switzerland,

RTCal<sup>™</sup> is a novel technology that provides heat flow profiles in real time.

RTCal<sup>™</sup> is complementary to the traditional heat flow method and adds additional data which translates into better information. The ease-of-use makes RTCal<sup>™</sup> a true walk-up tool that substantially enhances the capabilities and applications of the Reaction Calorimeter RC1e.

Measurements with RTCal<sup>TM</sup> are independent of the properties or the behavior of the reaction mass. Thus, investigation of previously difficult-to-study chemistries, such as polymerizations or reactions that may require immediate interaction, will become easier. The aspect of user-friendliness is further enhanced by iControl; the new reactor control and evaluation software which combines functionality with flexibility and a simple and straightforward graphical user interface.

# 1. Introduction

<u>RTCal<sup>TM</sup></u> is the new Real Time Calorimetry for the RC1e<sup>TM</sup> (Mettler-Toledo). The patented measurement technology <u>RTCal<sup>TM</sup></u> is based on heat flux sensors located in the jacket of a double-walled reactor vessel. Vertically and horizontally positioned sensor bands detect the heat flow Qrtc across the wall of the reactor online and in real time.

The heat flow information is sub-sequently processed, stored and directly displayed as Watts. Depending on the depth of information that you wish to obtain the system can additionally compute the overall heat of reaction Qr\_est including heats of dosing and accumulation and the integrated value IntQr\_est in real time.



*Figure 1*: *RTCal reactor with integrated sensors. The sensors consist of a horizontal band all around the reactor and a long sensor band in the back.* 

Experiments with RTCal<sup>TM</sup> are independent of the properties or the behavior of the reaction mass. Hence, no calibration procedures need to be run before, during or after the reaction which reduces the experiment time substantially compared with traditional methods.

The RTCal<sup>TM</sup> methodology can be run as a sole method on the RC1e<sup>TM</sup> (Mettler-Toledo) or complementary to the more traditional heat flow calorimetry (Visentin, F 2005). The enhanced set of information allows the user to derive more detailed conclusions and gives full confidence in experimental results.

The Ealphent   Calation Course (31, 2017(Analize)   PV	Contract Rice Contractory	
O iControl Rola-		METTLER TOLEDO
Experiment	Equipment	Perminders
New Experiment	Durige is view mis insument carigoration	The Experiment The Experiment Here a speciment
		양 Custod Holp 양 Custod in the wells 양 Product Catalog 양 License Activities Guide
Chemicals	Services	a magina
New Cherscal Manage Cherscals	Mixing Guidsfines Hardware Manuals	
Recard Documents		Contact the
NV Container, FTC at, IP No. 10 (11) (12) (12) (12) (12) (12) (12) (12)	Carry of Carrier Carry C	Sender-Lendtz-Centel Social Engl Control Control

*Figure 2: RTCal reactor in the new iControl RC1e software: Heat flow Qrtc, reaction heat Qr\_est, and its integration IntQr\_est are shown in real time during the experiment. Reactions can be optimized directly based on the information, e.g. by immediate changes of a dosing rate.* 

RTCal<sup>™</sup> is controlled and the results are evaluated by a new software, iControl (see Figure 1). The new graphic user interface improves the analysis in easy evaluation environment and shows all relevant data as trends or digital data at any time. The "Direct Mode" allows the user to immediate change all the parameter (stirring speed, temperature heat/cool rate, dosing, etc...) recording individual tasks and generate the procedure automatically (see Figure 3)



*Figure 3:* Screen-shoot of the experiment window showing the "Procedure", the "Equipment set-up" and the "Trend evaluation" during a running experiment.

## 2. Experimental section and results

The experiments were performed using a RC1e Reaction Calorimeter that combines two different calorimetric principles, heat flow and the new RTCal<sup>TM</sup>.

The sample volume of the reactor is 70 to 500 ml, and it is able to withstand pressure from 20 mmbar to 1 bar. The maximum temperature of the reactor is approximately 160  $^{\circ}$ C, and the minimum temperature depends only on the cryostat used down to -50 $^{\circ}$ C. Isothermal conditions are maintained using the RC1e internal thermostat (Riesen et. al. 1985, Mettler-Toledo).

The data acquisition is done using iControl software (see below).



*Figure 4*: Equipment set-up for the reactions: RC1e Reaction Calorimeter with a new 500 ml RTCal reactor run by the iControl RC1e software.

#### 2.1 Esterification reaction



Figure 5: Simplified reaction scheme for esterification reaction

Table 1: Recipe used for the esterification experiment.

Solvent:	Butanol		
Substrate:	Propionic anhydride		
Stirrer speed:	500 rpm <sup>a</sup>		
Temperature in the reactor:	70°C		

The standard measurements were made according to the following procedure. Firstly, the reactor was filled with butanol absolute and the stirrer was set at 500 rpm and the desired reaction temperature was set up to 70°C. In a second step the propionic anhydride was filled during 20 minutes.

To determine the energy introduced by dosing,  $q_{Dos}$ , the heat capacity of the feed (~2 kJ·kg<sup>-1</sup>·K<sup>-1</sup>) (NIST) was used. The reaction was followed for approximately three hours, until no further reaction was observed.



**Figure 6**: Comparison between the measured heat signal obtained with the standard method (heat flow) and the new RTCal<sup>TM</sup> method (Real Time Calorimetry). Reaction conditions: substrate: propionic anhydride; solvent: butanol; temperature:  $70^{\circ}C$ ; stirrer speed: 500 rpm.

Figure 6 shows the comparison between the heat flow measure with the standard method and the heat flow measured with the new RTCal<sup>TM</sup> principle. As can be seen the two trends are in good agreement.

The target of this esterification is the strong heat of dosing do to the addition of the propionic acid.

The reaction according to the conditions of Table 1 gives the following enthalpies. The integration of Qr could be done using the heat term given by the standard method  $Q_{flow}$  or using the new  $Q_{rtc}$  signal. The numerical results are shown in Table 2.

Experiment integration based on	Qr	Qflow	Qrtc	Deviation (%)
Esterification integration results	-67.4 kJ/mol	-54.9 kJ/mol	-56.9 kJ/mol	+ 3.5

Table 2: Reaction enthalpies for the esterification experiment.

This high consistency demonstrates the high accuracy of the measurements that come from the two different methods.

2.2 Grignard Reaction



Figure 7: Simplified reaction scheme for the Grignard experiment.

Table 3: Recipe used for the Grignard reaction

Solvent:	THF + toluene
Reagent 1:	Magnesium
Reagent 2:	Parabromotoluene + toluene
Stirrer speed:	500 rpm <sup>a</sup>
Temperature in the reactor:	70°C

The standard measurements were made according to the following procedure. Firstly, the reactor was filled with THF, toluene and magnesium and the stirrer was set at 500 rpm and the desired reaction temperature was set up to 40°C. In a second step the parabromotoluene and toluene was filled during 20 minutes.

To determine the energy introduced by dosing,  $q_{Dos}$ , the heat capacity of the feed (~2 kJ·kg<sup>-1</sup>·K<sup>-1</sup>) (NIST) was used. The reaction was followed for approximately three hours, until no further reaction was observed.

The reaction was started with THF, Toluene, and Magnesium (118g in total) at  $40^{\circ}$ C. The temperature was kept constant by dosing Parabromtoluene and Toluene (196 g in total) into the reactor. The resulting reaction is shown below:



*Figure 8*: grignard reaction monitored in real time: The blue curve shows the mass of the reaction mixture mr, i.e. the dosing of parabromtoluene and toluene. In green the heat of reaction obtained in real time based on RTCalTM, in red the reaction temperature Tr.

As Figure 8 shows,  $RTCal^{TM}$  monitors a very strong peak of reaction heat at the beginning of the reaction up to Qr = 150W in the reaction mixture. The reaction enthalpy resulting in real time fits nicely with former values of the standard method as shown in the table below.

Table 4: Reaction enthalpies for the Grignard experiment.

Experiment integration based on	Qr	Qrtc
Grignard integration results	-510 kJ/mol	-487 kJ/mol

#### **3.** Conclusions

The new RTCal<sup>TM</sup> principle has been used to characterize two reactions examples. The esterification of propionic anhydride was performed in order to demonstrate the feasibility, precision and advantage of following the reaction with a strong heat of dosing. The Grignard reaction demonstrate instead that RTCal<sup>TM</sup> could be used for safety studies to characterize and optimize in real time dangerous reactions.

RTCal<sup>TM</sup> heat flow profiles therefore allows the user to make adjustments to the actual chemistry and optimization of process parameters online.

RTCal<sup>TM</sup> is complementary to the traditional heat flow method and adds additional data which translates into better information. The ease-of-use makes RTCal<sup>TM</sup> a true walk-up tool that substantially enhances the capabilities and applications of the Reaction Calorimeter RC1e<sup>TM</sup> (Mettler-Toledo).

iControl is the new evaluation and control software which combines functionality with flexibility and a simple and straightforward graphical user interface. This modern and state-of the-art user interface reflects the personal way to run experiments and offers what was always expected from the control software.

In summary, regardless of whether you are looking for immediate thermal data, searching for specific details within your process, generating data to develop a kinetic model or simply optimizing a process RTCal<sup>TM</sup> will help to find the answer in an efficient and cost effective manner.

### 4. References

- Riesen R., Grob B., Reaktionskalorimetrie in der chemischenProzessentwicklung, Swiss Chem, Vol. 7, No. 5a, pp. 39-43, 1985
- Mettler-Toledo, www.mt.com
- NIST Chemistry Webbook, Standard Reference Database Number 69, National Institute of Standards and Technology, online, 2003
- Visentin, F. Ph.D. Thesis, Kinetic Study of Hydrogenation Reactions of Aromatic Nitro Compounds Using a New Pressure Resistant Reaction Calorimeter Combined with a FTIR-ATR Device. In Diss. ETH No. 16053, Zurich, Switzerland, 2005.